


Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

XRF Instrumentation

Introduction to sources


AMPTEK, INC.

14 Deangelo Drive, Bedford, MA 01730
Ph: +1 781 275 2242 Fax: +1 781 275 3470
sales@amptek.com www.amptek.com


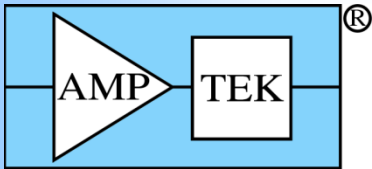


Instrument

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

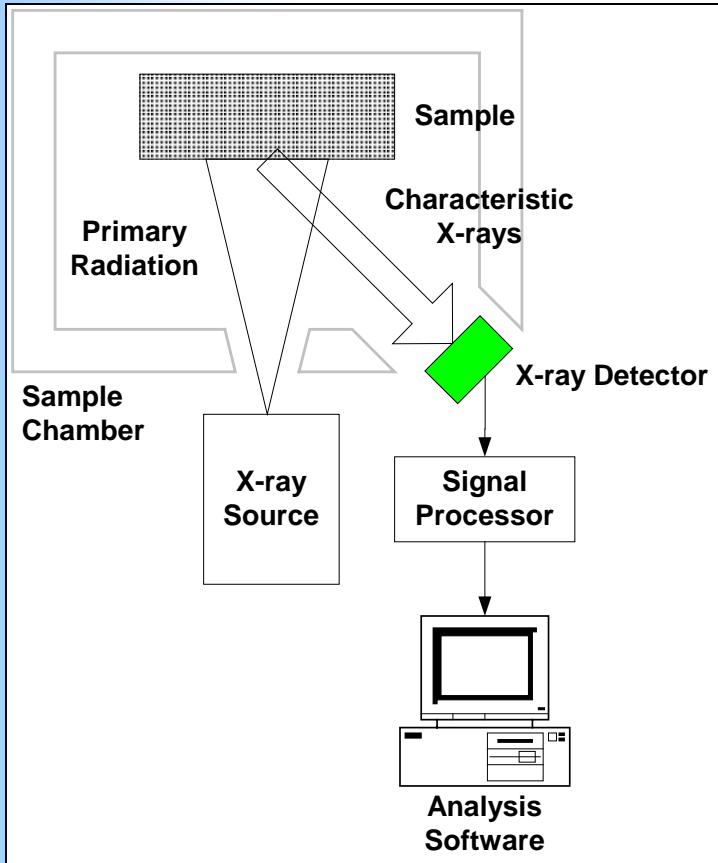


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- Excitation source
 - X-ray tube or radioisotope
- Spectrometer
 - X-ray detector
 - Signal processing electronics
- Software
 - Spectrum correction and processing software
- Other
 - Radiation shielding
 - Sample fixture

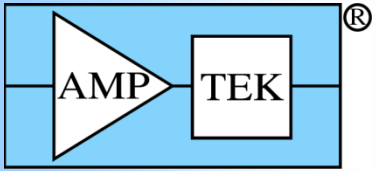

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Excitation Source

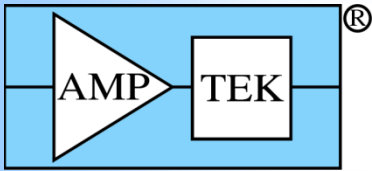

Sources



Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

What is the purpose of the source?

- Excite elements to be analyzed
- Produces the signal


What is an ideal XRF source?


- Energy just above edge of analyte
- Monoenergetic → Low background
- Intense flux → Fast measurement
- Stable flux → Accuracy
- Can be turned off → Radiation safety
- Long life, inexpensive

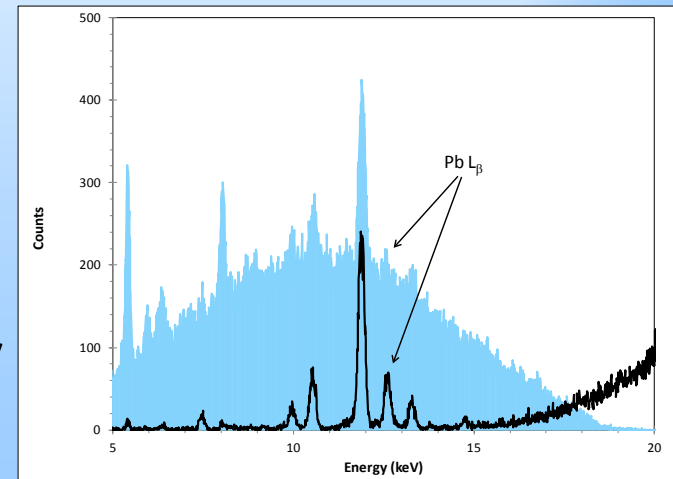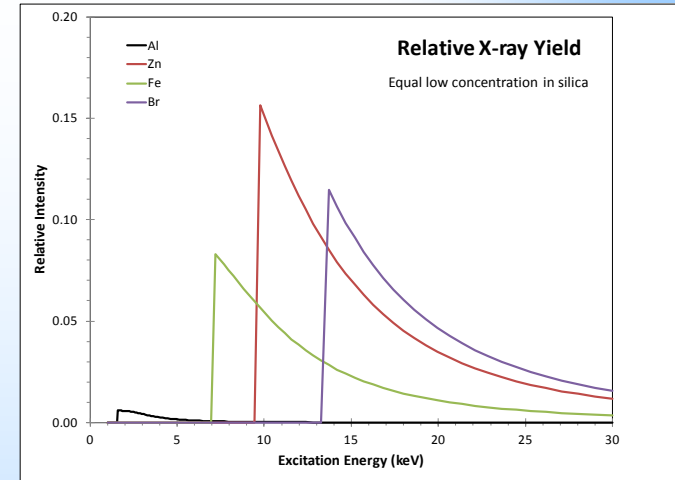

X-ray tubes

- Heated filament produces electrons
- High voltage (10 to 50 kV) target accelerates electrons
- Electrons hitting target → X-rays
 - Characteristic X-ray lines from target material
 - Bremsstrahlung continuum from electron scattering
- Many variations
 - Anode material
 - Maximum bias and current
 - Side window, end window, transmission
 - Microfocus, pulsed, DC

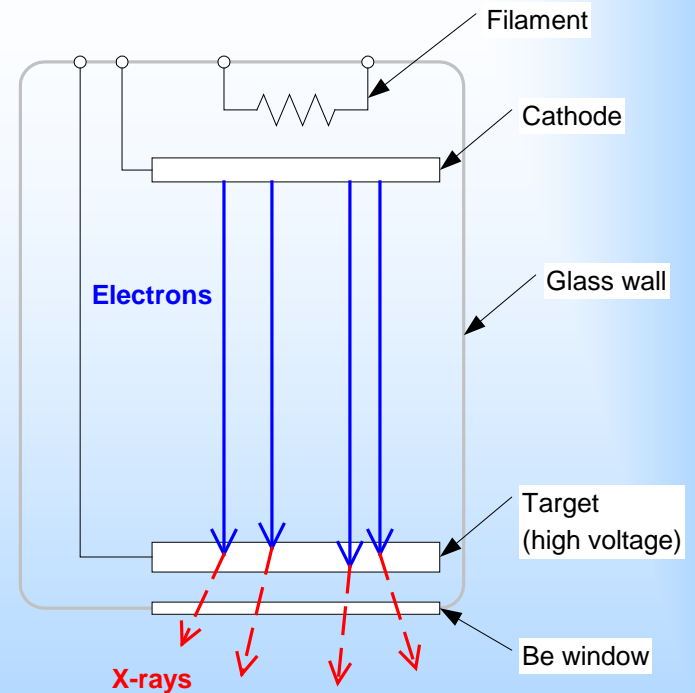

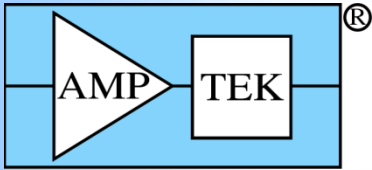


Sources

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

X-ray tube variations

- Geometry
 - End window: Permits a very tight geometry but power is limited
 - Side window: Allows much higher power but also much larger
- Anode material
 - Produces characteristic X-rays. Can be useful for exciting sample or can interfere
 - Ag is very common (almost a standard). Rh, Pd, Au, W also used
- Maximum HV and current
 - HV setting is critical for exciting sample. High current permits fast measurement
 - Higher HV and current require more shielding and radiation safety issues
- Specialties
 - Microfocus
 - Pulsed

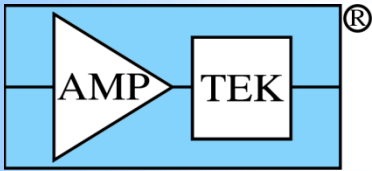



Sources

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

X-ray tubes for XRF

- Disadvantages
 - Most X-rays are in continuum → High background
 - Large, high power, heat dissipation
- Advantages
 - Can change HV and filters → Tune excitation for different elements
 - High intensity possible
 - Can turn it off, improves safety

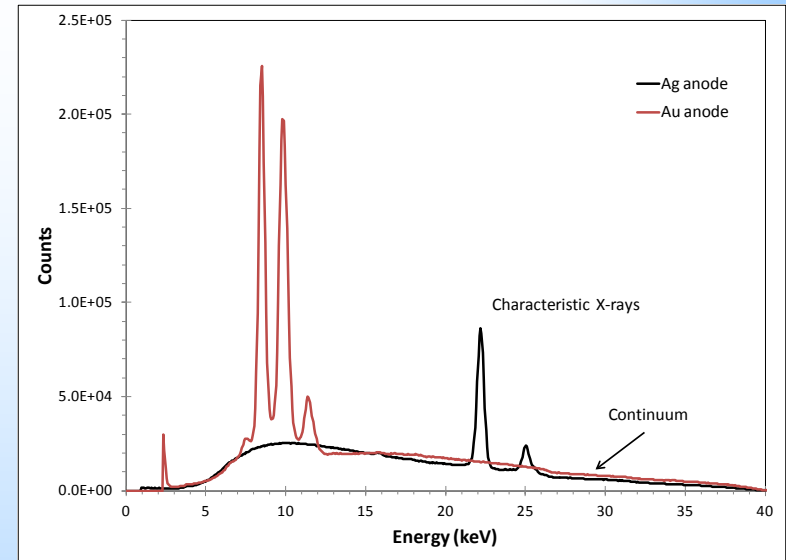

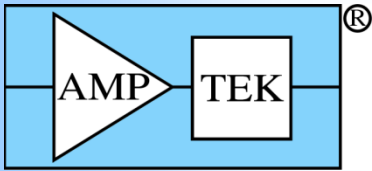

®

Sources


Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Radioisotopes

- Monoenergetic → Low background
- No power dissipation
- Fixed, limited emission energies
- Cannot turn them off
- Intensity falls with half life

- Modern tube technology has replaced sources in many but not all applications

	^{55}Fe	^{109}Cd	^{241}Am	^{57}Co
Energy (keV)	5.9	22.1 88	59.5	122
Elements (K lines)	Al - V	Fe - Mo	Ru - Er	Ba - U
Elements (L lines)	Br - I	Yb - Pu		



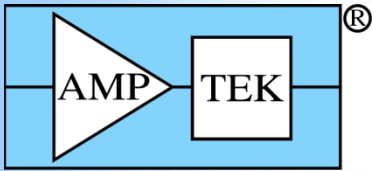



Sources

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Source filter

- Attenuates incident X-ray below attenuation edge of filter material
- Reduces background
- Greatly enhances signal to background
- Optimum filter depends on energy of X-rays to be analyzed

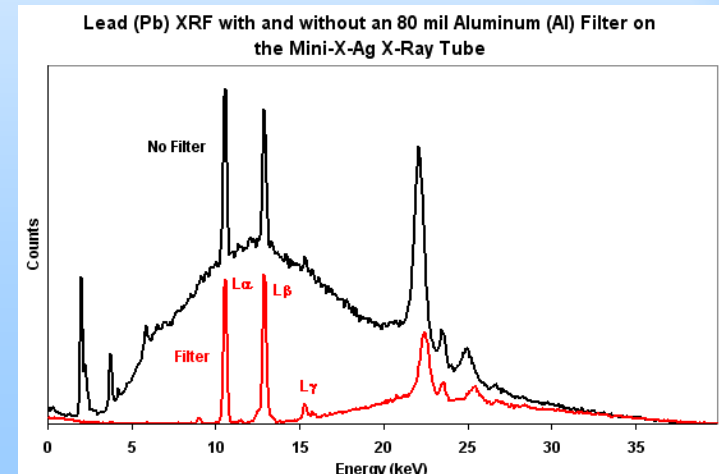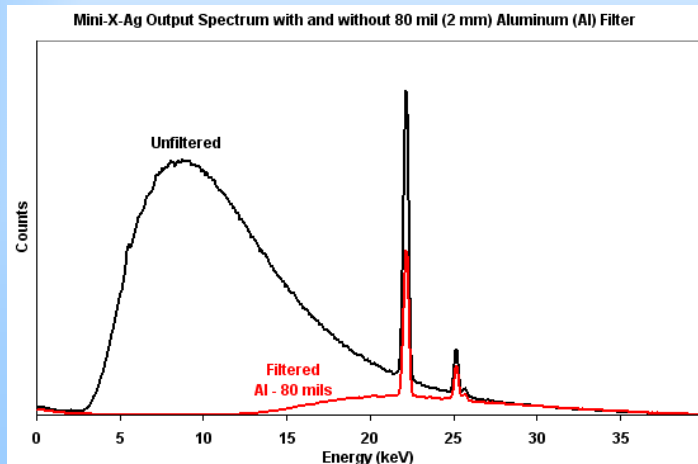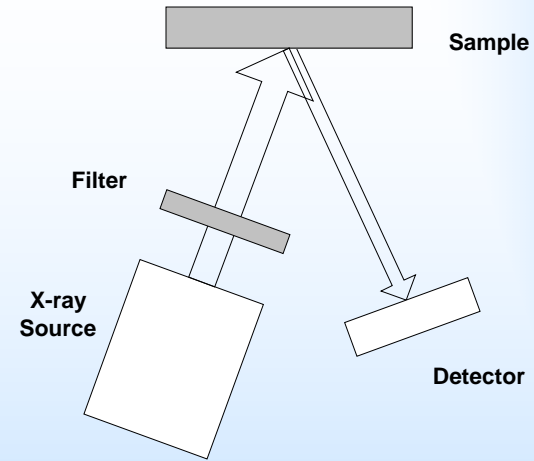

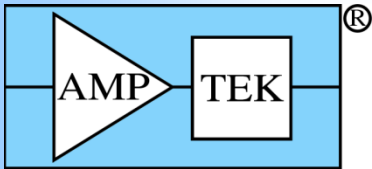


Sources - Advanced

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Secondary targets

- Primary beam excites K lines of the secondary target, which excite the sample.
- Excitation is at target K lines, so background is greatly reduced
- Inefficiency of the process requires much higher X-ray tube power.

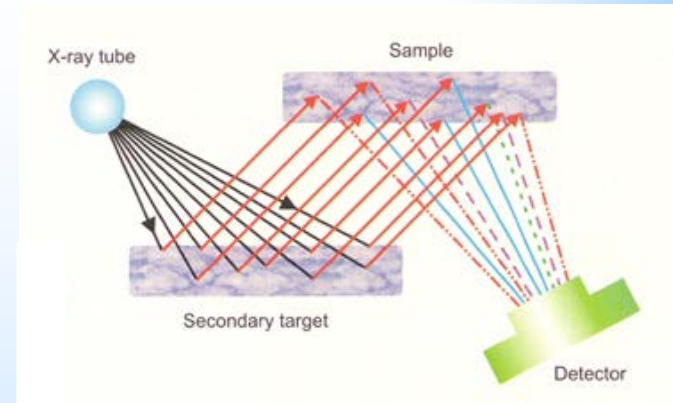



Capillary Optics

- X-rays cannot be reflected or refracted but capillary optics can be used to concentrate
- Increases signal intensity

Crystal Optics

- Crystals are used to diffract primary beam
- Only transmits in a narrow energy range so reduces background

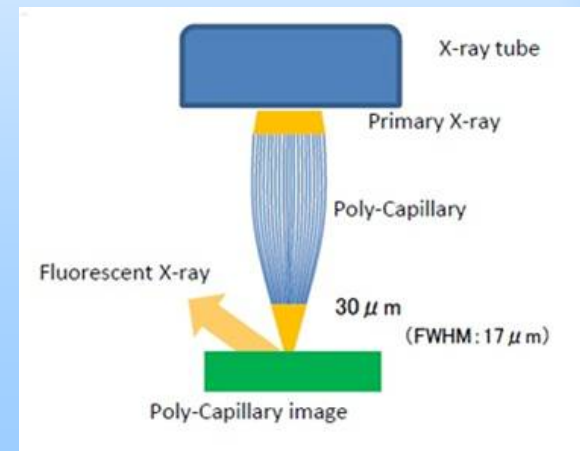

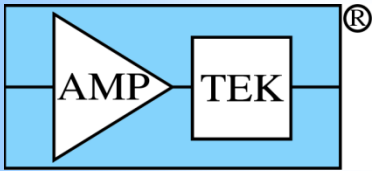

Sources - Advanced

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Electron beams


- Also excite characteristic X-rays. Used in "electron dispersive spectroscopy", related to XRF but different method
- Main advantage is that electrons have short range in sample so are very effective for measuring low energy X-rays from the lightest elements

Particle induced X-ray emissions (PIXE)

- Alpha particles, protons, and other energetic particles ionize materials which thus give off characteristic X-rays
- Can be used for X-ray spectroscopy (though not strictly XRF)

Synchrotrons

- Extremely bright source of X-rays so permit sophisticated analysis
- Only available at these large research facilities

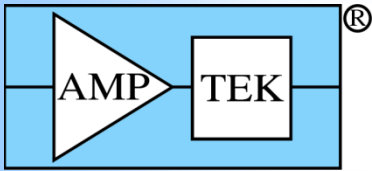


Amptek X-ray Sources

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Mini-X and Mini-X OEM

- 50 kVp, 4W, small tube with USB control
- Au, Ag & Rh anodes, 10W tubes available
- Easy to attach custom collimator



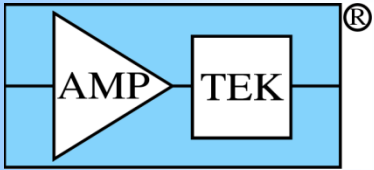

AmpTek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Additional Advice


Novice users often underestimate the important of optimizing the excitation source. They assume that, if they are exciting X-rays, the spectrometer is configured properly, and the analysis software has been set right, they will get good results.

One will, indeed, be able to see a signal. But the precision, accuracy, and detection limits will be much worse than they anticipate unless they are careful with the excitation conditions.

Rules of thumb:

- Excitation energy must be 2 keV above line you are measuring to see anything
- Best excitation is usually at a kV 2-3 times the energy of the line
- Best filter will use a K edge about 1.5 time energy of the line

40 kVp with a 10 mil Al filter is "jack of all trades, master of none". It will show all lines but will measure none of them very well.


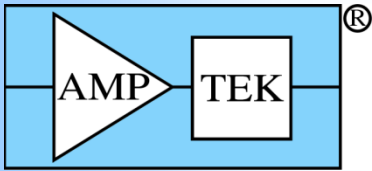



Sources

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Filters and kVp are key

- Plots below show same targets with different filters
- Blue: No filter, good signal at low energy
- Red: Filter gives much better signal to background

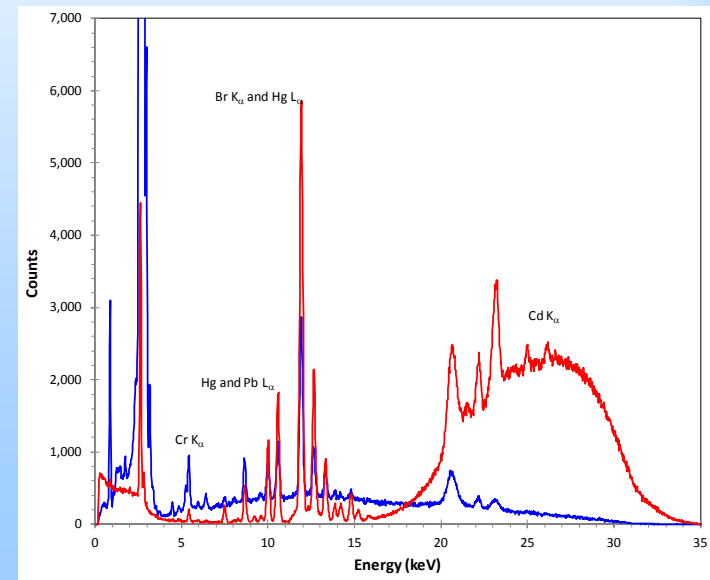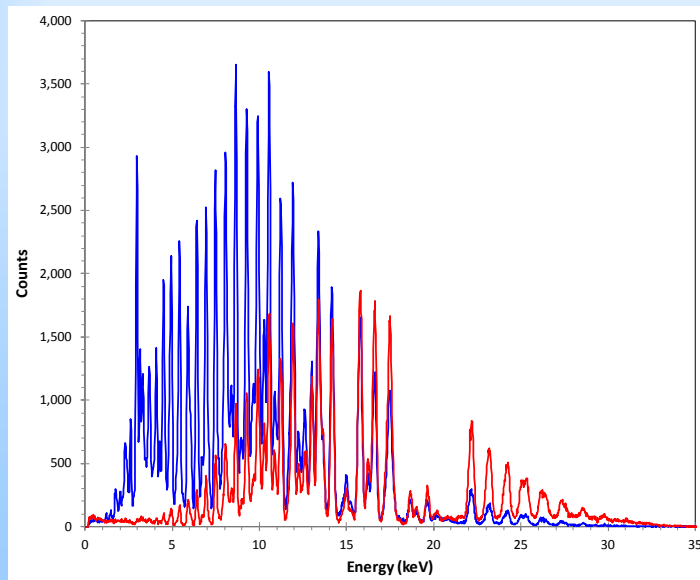


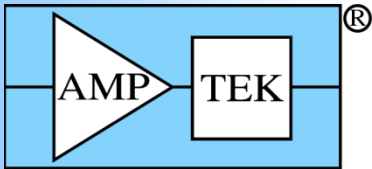



Sources

Amptek Inc.
14 Deangelo Drive
Bedford, MA 01730 USA
www.amptek.com

Filters and kVp are key

- Plots here show three options
- W/Al filter reduces 5 to 10 keV background x500
- X-rays penetrate deeply so little signal < 5 keV
- No filter only best at the lowest energies

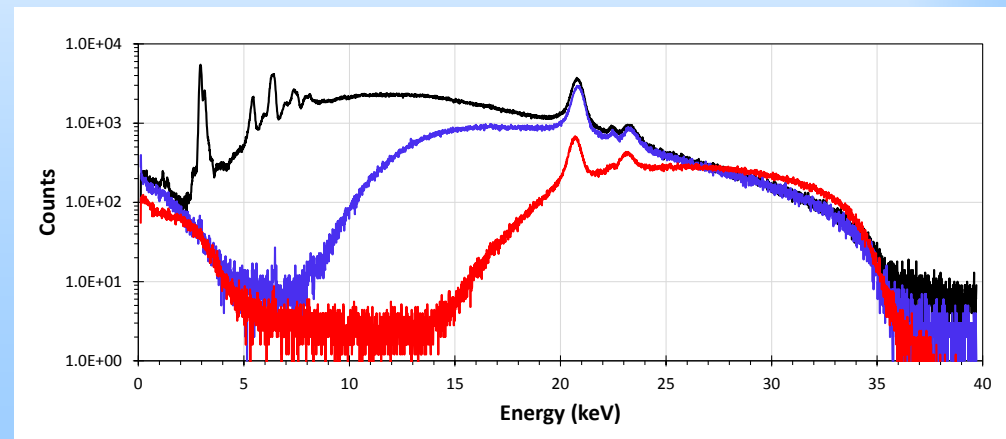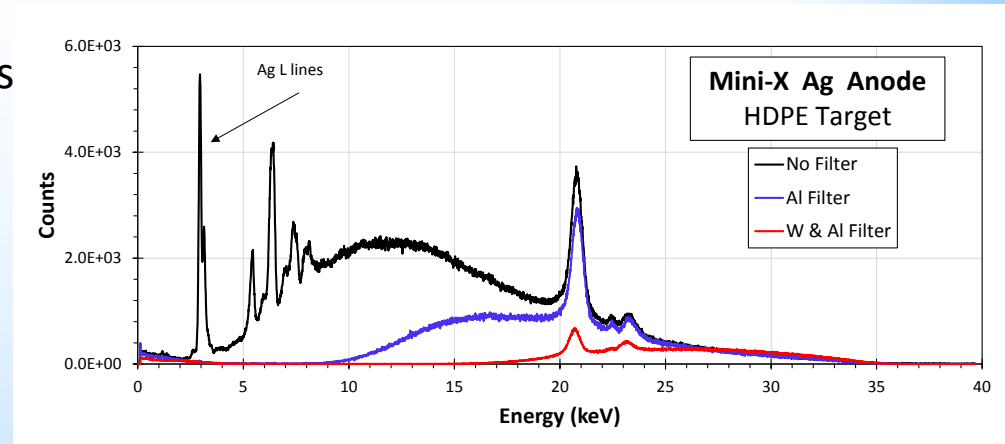

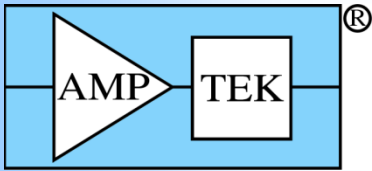

Sources

Amptek Inc.
 14 Deangelo Drive
 Bedford, MA 01730 USA
 www.amptek.com

Filters and kVp are key

- Different filter and kVp setting gave best results for a different range of elements

	No Filter 15 kVp		No Filter 25 kVp		No Filter 35 kVp		No Filter 45 kVp		Filter 35 kVp		Filter 45 kVp	
	σ	S/B	σ	S/B	σ	S/B	σ	S/B	σ	S/B	σ	S/B
Ti	2.05%	2.80	2.44%	2.94	2.86%	2.77	2.82%	3.08	15.47%	1.56	21.71%	0.78
Fe	1.55%	5.48	1.68%	6.21	1.86%	6.31	1.98%	6.05	5.62%	6.24	6.18%	4.61
Ni	1.63%	4.97	1.58%	6.02	1.65%	6.69	1.76%	6.42	3.51%	8.32	3.68%	7.47
Zn	2.50%	1.98	8.12%	0.46	2.78%	1.73	5.11%	0.86	5.13%	1.97	4.47%	2.33
Ga	2.35%	2.16	1.73%	2.70	1.73%	2.79	1.82%	2.64	2.95%	2.91	3.02%	2.79
Ge	2.88%	1.92	1.83%	2.53	1.76%	2.61	1.78%	2.59	2.76%	2.65	2.81%	2.62
As	17.81%	0.28	1.86%	2.64	1.75%	2.76	1.73%	2.81	2.40%	3.17	2.40%	3.18
Se	6.93%	1.09	2.24%	3.27	2.87%	2.23	1.95%	3.84	2.79%	4.30	3.20%	3.39
Br	24.28%	0.30	1.93%	3.96	1.92%	3.59	1.72%	4.01	2.14%	4.76	2.31%	4.10
Rb			2.51%	3.19	1.99%	3.61	1.77%	4.06	1.90%	4.98	2.03%	4.36
Sr			2.35%	5.64	1.89%	6.94	1.52%	9.95	1.67%	13.53	1.97%	10.07
Zr			3.40%	3.21	2.44%	4.07	1.70%	6.42	1.64%	8.56	2.11%	5.58
Mo			4.12%	3.61	2.12%	7.78	1.65%	11.29	1.43%	19.42	1.66%	14.41
Ag					5.29%	3.13	4.16%	4.28	2.32%	6.16	2.52%	6.27
Sn					17.95%	0.81	4.40%	4.55	7.80%	0.91	4.17%	2.19